Fig. 1

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

## UK – System 50 TM

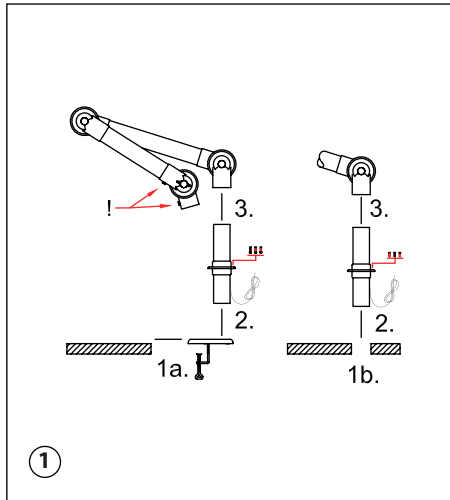**Fig. 1**

On account of the packaging the extraction arm is delivered in two parts.

1. For mounting in an ESD working area the earthing wire on the socket pipe is very important. Take care not to damage it during the mounting.
2. If the table is of a conductive material the connection to the extraction arm/bracket must be non-conductive.
3. Mount the table bracket on the edge of the table and tighten the clamps (fig. 1a). Or make a hole in the table with a diameter of 60mm where the extraction arm is to be placed (fig. 1b).
4. Place the socket pipe in the hole of the bracket/table and fasten the flange to the bracket/table (2).
5. Place the joint of the extraction arm on the socket pipe. Be aware that the spring goes into the pipe (3) and the finger-screws (!) face forward.

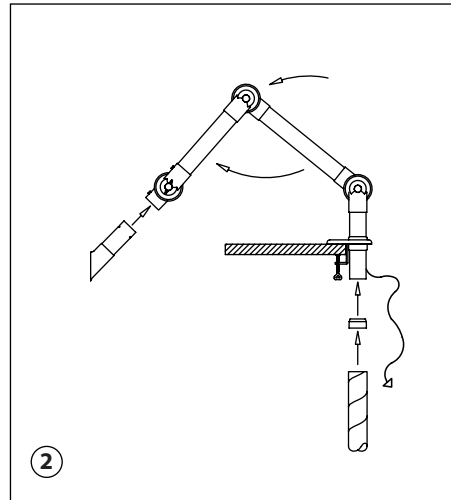
**Fig. 2**

1. Connect the extraction arm to the main ventilation duct with the Alsident® reducer.
2. In an ESD area a potential equalizing must be ensured for the surrounding working area. The earthing wire with the 1MΩ resistance must be connected to the central earthing for the furniture and tools. Take care not to break the 1MΩ resistance.
3. Unfold the extraction arm in the direction of the arrows.
4. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small fingerscrew on the joint. When screwed in the fingerscrew will hold the accessory in place in the joint.
5. Finally make sure the damper is open.
6. The extraction arm is now ready for use.

Alsident® System reserve the right to make changes in this brochure to technical data, construction, materials, equipment and appearance. The illustrations do not necessarily represent the true appearance of the products and are for the purpose of description only. Alsident® System make reservations regarding misprints.

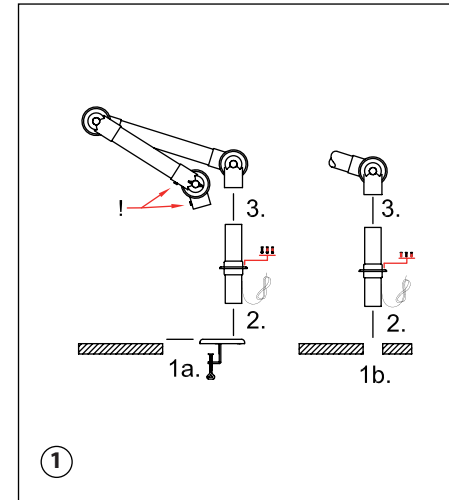See MAINTENANCE AND CLEANING at [www.alsident.com](http://www.alsident.com) under "Technical Information"

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