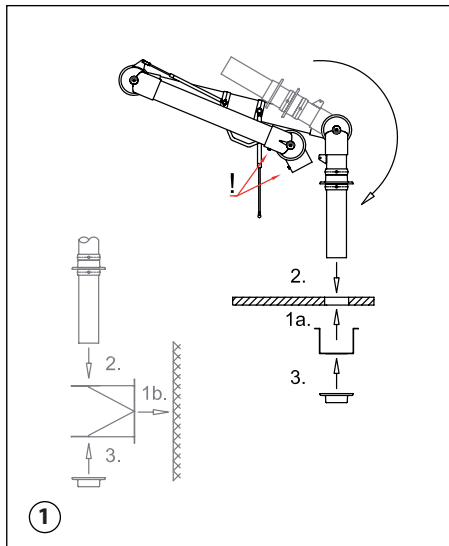
## Alsident® System 100

Aluminium and Chemical Resistant  
Aluminium und Chemisch Beständig  
Aluminium og Kemisk Resistent

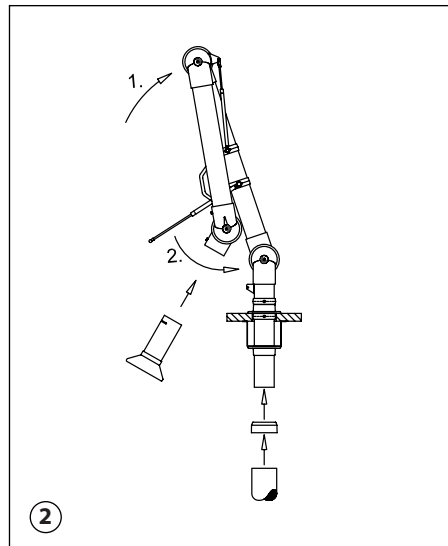
**Fig. 1**

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
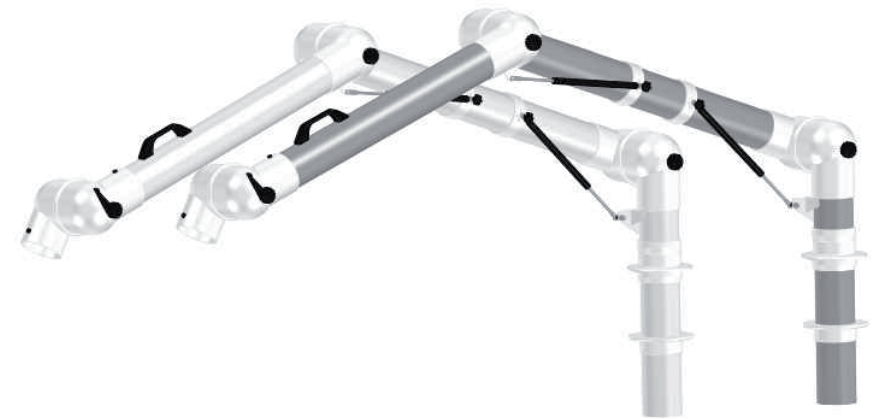
## UK – System 100 TM Gas spring

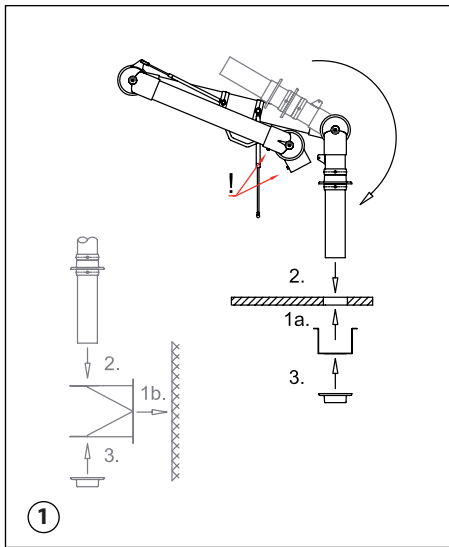
**Fig. 1**

1. Make a hole with a diameter of 120 mm where the extraction arm is to be placed. Fasten the U-profile to the table (1a). Or fasten the wall bracket on the wall in the desired height above the floor (1b).
2. When unfolding the extraction arm, turn the socket pipe in the direction of the arrow. Make sure the fingerscrews (!) are faces forward.
3. Place the socket pipe through the holes in the U-profile/wall bracket and fasten the flange with screws (2).
4. Place the last flange on the socket pipe under the table/bracket and fasten it with screws (3).

*NB! Gas springs are considered as wearing parts.*

Alsident® System reserve the right to make changes in this brochure to technical data, construction, materials, equipment and appearance. The illustrations do not necessarily represent the true appearance of the products and are for the purpose of description only. Alsident® System make reservations regarding misprints.

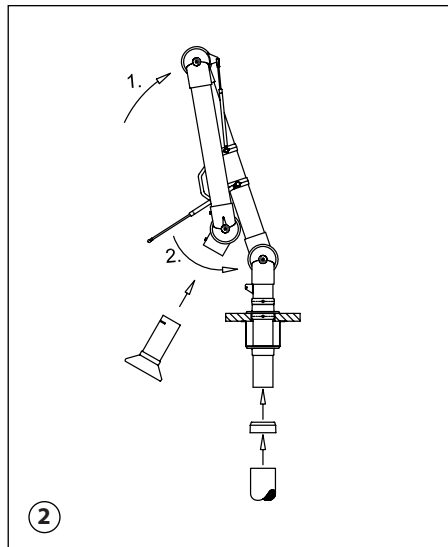
**Fig. 2**

1. To mount the gas spring turn the extraction arm towards vertical position (1)
2. Turn the gas spring to the fitting on the socket pipe. Fasten the gas spring to the fitting (2).
3. Connect the extraction arm to the main ventilation duct with the Alsident® reducer. In order to turn the extraction arm, it is important not to fix the reducer to the socket pipe.
4. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small fingerscrew. When screwed in the fingerscrew will hold the accessory in place in the joint.
5. Finally make sure the damper is open.
6. The extraction arm is now ready for use.

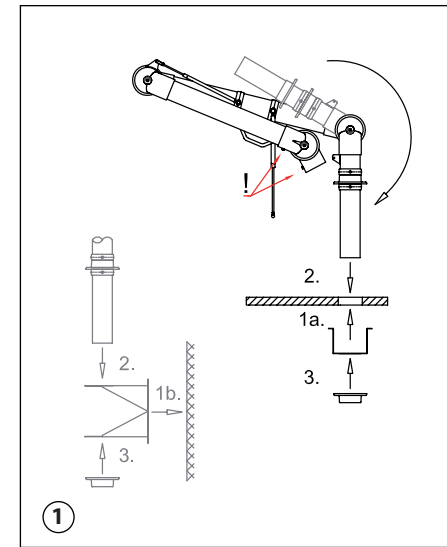See MAINTENANCE AND CLEANING at [www.alsident.com](http://www.alsident.com) under "Technical Information"

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