BM  
gas

# Alsident® System 100

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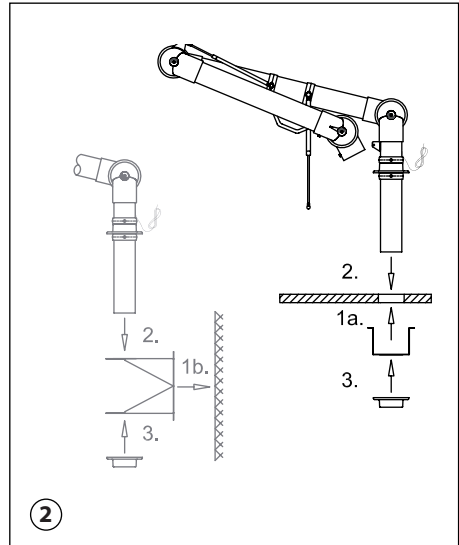
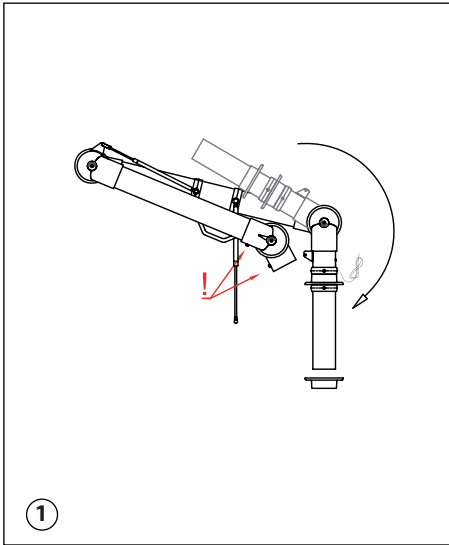
UK

**Fig. 1 (EX/ESD):**

1. When unfolding the extraction arm the socket pipe is turned in the direction of the arrow. Make sure the fingerscrews (!) are faces forward.
2. For mounting in an ESD/EX working area the earthing wire on the socket pipe is very important. Take care not to damage it during the mounting.

**Fig. 2 (EX/ESD):**

Please note! If the table is of conductive material the connection must be non-conductive.

- 1a. Make a hole with a diameter of 120 mm where the extraction arm is to be placed. The U-profile is fastened to the table (above or below).
- 1b. The wall bracket is fastened on the wall or column in the desired height above floor.
2. The socket pipe is placed through the holes in the U-profile/wall bracket and the flange is fastened with screws.
3. Place the last flange on the socket pipe under the table/bracket and fasten it with screws.

*NB! Gas springs are considered as wearing parts.*

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
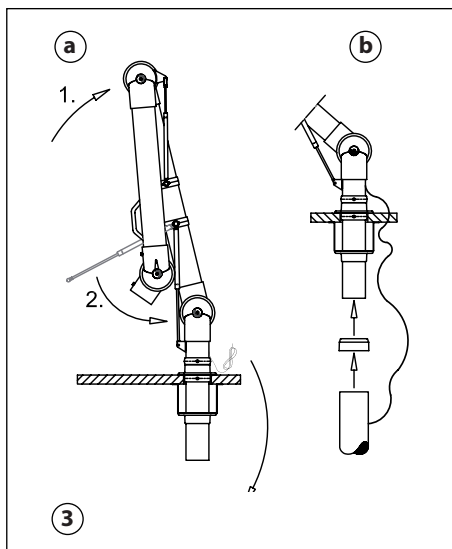
UK

**Fig. 3 (EX):**

1. To mount the gas spring turn the extraction arm towards vertical position. (fig. 3a)
2. Turn the gas spring to the fitting on the socket pipe. Fasten the gas spring to the fitting.
3. Connect the extraction arm to the main ventilation duct with the Alsident® reducer (fig. 3b).
4. In an explosive working area a potential equalizing must be ensured for the ventilation system. It is therefore important to bond all the conductive parts of the system and earthing them. The extraction arm must be connected to the ventilation duct with the earthing wire pre-attached to the extraction arm. (fig. 3.b). Alsident® System A/S recommends the earthing of the extraction arm to be carried out by personnel experienced in EX areas.

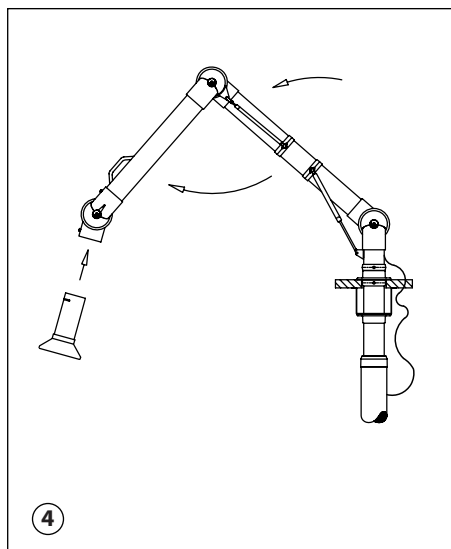**Fig. 4 (EX):**

1. Unfold the extraction arm in the direction of the arrows.
2. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small fingerscrew. When screwed in the fingerscrew will hold the accessory in place in the joint.
3. Finally make sure the damper is open.
4. The extraction arm is now ready for use.

*Approved for use in explosive areas and marked*

**EX II 1 GD**

*according to standards: IEC 61340-4-1:2003 and  
EN 13463-1:2001, supporting ATEX directive 94/9/EC.  
For further information see Technical Information at:  
[www.alsident.com](http://www.alsident.com).*

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
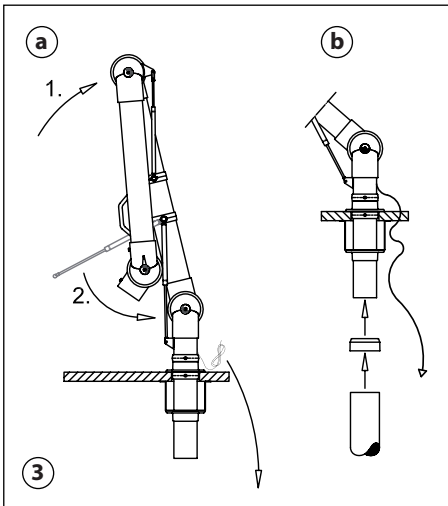
UK

**Fig. 3 (ESD):**

1. To mount the gas spring turn the extraction arm towards vertical position (fig. 3a).
2. Turn the gas spring to the fitting on the socket pipe. Fasten the gas spring to the fitting.
3. Connect the extraction arm to the main ventilation duct with the Alsident® reducer (fig. 3b).
4. In an ESD area a potential equalizing must be ensured for the surrounding working area. The earthing wire pre-attached to the extraction arm must be replaced with the one in the plastic bag marked "ESD". The new earthing wire with a 1MΩ resistance must be connected to the central earthing for the furniture and tools (fig. 3b). When turning the extraction arm be aware not to break the 1MΩ resistance on the earthing wire. Be aware to break any conductive contact between the bracket and the building/furniture and the main ventilation system.

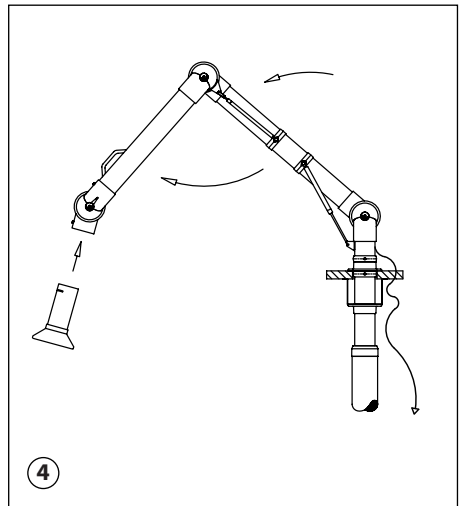**Fig. 4 (ESD):**

1. Unfold the extraction arm in the direction of the arrows.
2. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small fingerscrew. When screwed in the fingerscrew will hold the accessory in place in the joint.
3. Finally make sure the damper is open.
4. The extraction arm is now ready for use.

Approved according to standard IEC 61340-5-1:2003.  
For further information, see Technical Information at:  
[www.alsident.com](http://www.alsident.com).

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UK:

Alsident® System reserve the right to make changes in this brochure to technical data, construction, materials, equipment and appearance. The illustrations do not necessarily represent the true appearance of the products and are for the purpose of description only. Alsident® System make reservations regarding misprints.

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