HMS
gas

Alsident® System 100

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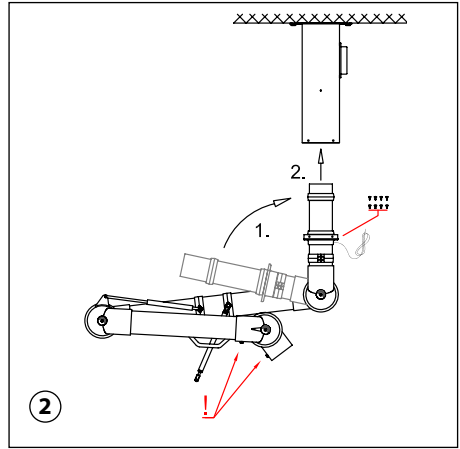
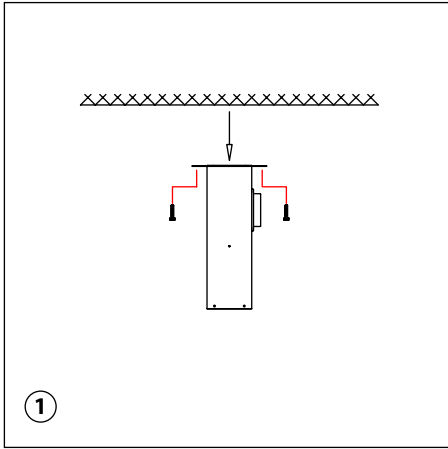
UK

Fig. 1 (EX/ESD):

1. Mount the column solidly on the ceiling. Be aware of the considerable strain the operation of the arm will transmit to the column.
2. The operation of the extraction arm will transmit a considerable strain to the column, therefore it must be mounted on a solid ceiling.

Fig. 2 (EX/ESD):

For mounting in an ESD/EX working area the earthing wire on the socket pipe is very important. Take care not to damage it during the mounting.

1. When unfolding the extraction arm, make sure the socket pipe is turned in the direction of the arrow and the finger-screws (!) are faced forward.
2. Place the socket pipe in the column. Turn the guide on the socket pipe until it falls into place in the column and fasten with the included screws.
3. The extraction arm can now be turned in the column.

NB! Gas springs are considered as wearing parts.

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
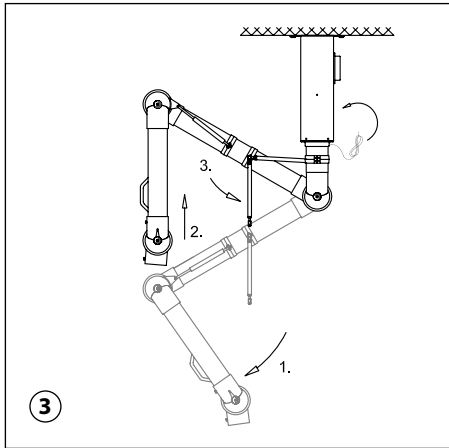
UK

Fig. 3 (EX):

1. Unfold the extraction arm in the direction of the arrow (1).
2. Push the extraction arm towards vertical position. This enables the gas spring to be mounted (2).
3. Turn the gas spring up towards the fitting on the socket pipe and fasten it with the finger-screw (3).
4. In an explosive working area a potential equalizing must be ensured for the ventilation system. It is therefore important to bond all the conductive parts of the system and earthing them. The extraction arm must be connected to the ceiling column with the earthing wire pre-attached to the extraction arm. The installer of the ventilation duct must make sure to connect the ceiling column to the ventilation duct with an earthing wire. (fig.4a).

Alsident® System A/S recommends the earthing of the extraction arm to be carried out by personnel experienced in EX areas.

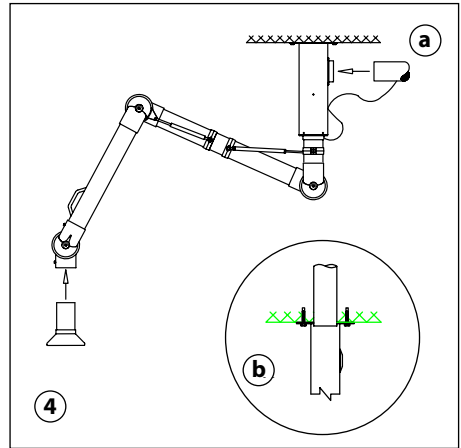**Fig. 4 (EX):**

1. The main ventilation duct (100 mm dia.) can be connected directly to outlet on the side of the column. (fig.4a). For top connection the main ventilation duct is put through the ceiling and directly into the column. Make sure to tighten the gap between the ceiling and duct with a filler. (fig. 4b).
2. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small finger-screw. When screwed in the fingerscrew will hold the accessory in place in the joint.
3. Finally make sure the damper is open.
4. The extraction arm is now ready for use.

Approved for use in explosive areas and marked

EX II 1 GD

according to standards: IEC 61340-4-1:2003 and EN 13463-1:2001, supporting ATEX directive 94/9/EC. For further information see Technical Information at: www.alsident.com.

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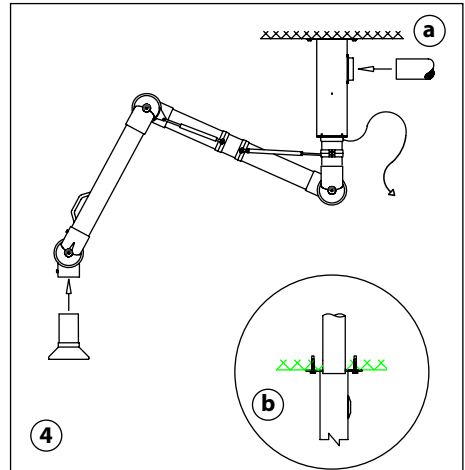
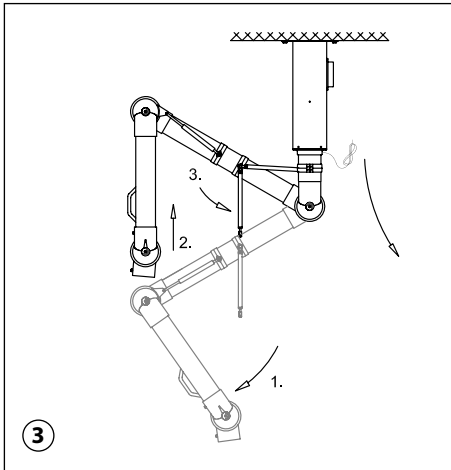
UK

Fig. 3 (ESD):

1. Unfold the extraction arm in the direction of the arrow (1).
2. Push the extraction arm towards vertical position. This enables the gas spring to be mounted (2).
3. Turn the gas spring up towards the fitting on the socket pipe and fasten it with the fingerscrew (3).
4. In an ESD area a potential equalizing must be ensured for the surrounding working area. The earthing wire pre-attached to the extraction arm must be replaced with the one in the plastic bag marked »ESD«. The new earthing wire with a 1MΩ resistance must be connected to the central earthing for the furniture and tools (fig. 4). When turning the extraction arm be aware not to break the 1MΩ resistance on the earthing wire. Be aware that there is no conductive connection to the main ventilations system.

Fig. 4 (ESD):

1. The main ventilation duct (100 mm dia.) can be connected directly to outlet on the side of the column. (fig. 4a).
For top connection the main ventilation duct is put through the ceiling and directly into the column. Make sure to tighten the gap between the ceiling and duct with a filler. (fig. 4b).
2. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small finger-screw. When screwed in the fingerscrew will hold the accessory in place in the joint.
3. Finally make sure the damper is open.
4. The extraction arm is now ready for use.

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UK:

Alsident® System reserve the right to make changes in this brochure to technical data, construction, materials, equipment and appearance. The illustrations do not necessarily represent the true appearance of the products and are for the purpose of description only. Alsident® System make reservations regarding misprints.

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