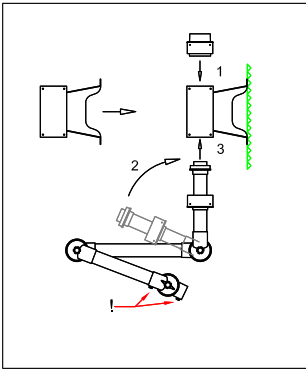
# WALL MOUNTED – WANDMONTAGE – VÆGMONTAGE


## UK – System 50 WM

**Fig. 1**

For mounting in an ESD working area the earthing wire on the socket pipe is very important. Take care not to damage it during the mounting.


1. Fasten the wall bracket solidly to the wall or column. When mounting the wall bracket be aware of the considerable strain the operation of the arm will transmit to the bracket.
2. If the wall or column is made of a conductive material the connection to the wall bracket must be non-conductive.
3. Place square-guide in the top of the bracket (1) and fasten it to the wall bracket with the included screws.
4. Considering the internal spring, it is important to turn the socket pipe in the direction of the arrow (2) and that the fingerscrews (1) face forward.
5. Place the socket pipe in the wall bracket (3) and fasten the square-guide with the included screws.
6. The extraction arm can be turned 190° in the wall bracket.

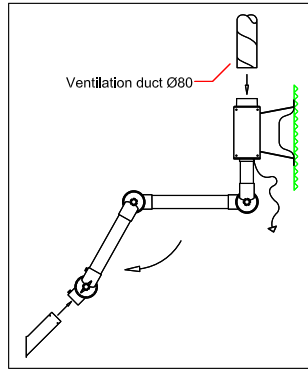**Fig. 2**

1. Connect the main ventilation duct (80 mm dia.) directly to the top of the wall bracket.
2. In an ESD area a potential equalizing work area. The earthing wire with the 1MΩ resistance must be connected to the central earthing for the furniture and tools. Take care not to break the 1MΩ resistance.
3. Unfold the extraction arm in the direction of the arrow.
4. Place the connection pipe of the accessory in the joint. Turn the slit in the connection pipe towards the small fingerscrew. When screwed in the fingerscrew will hold the accessory in place in the joint.
5. Finally make sure the damper is open.
6. The extraction arm is now ready for use.

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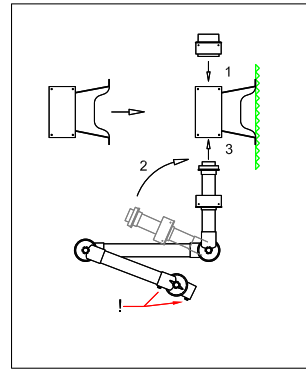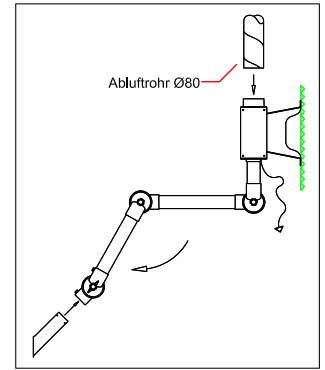
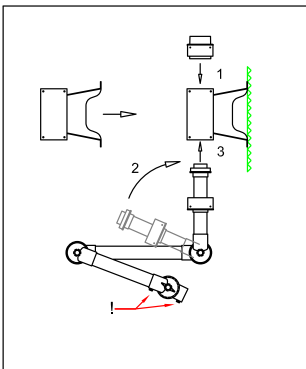
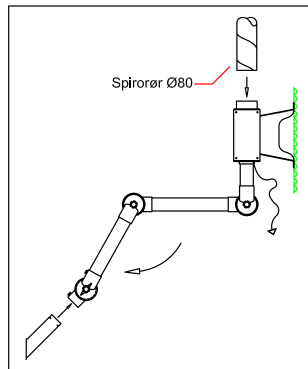
See MAINTENANCE AND CLEANING  
at [www.alsident.com](http://www.alsident.com) under "Technical Information"

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

## UK – System 50 CM

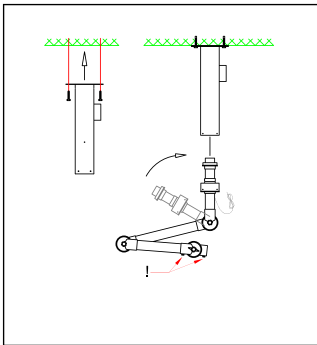**Fig. 1**

For mounting in an ESD working area the earthing wire on the socket pipe is very important. Take care not to damage it during the mounting.


1. Mount the column directly under the ceiling.
2. The operation of the extraction arm will transmit a considerable strain to the column. The column must therefore be mounted on a solid ceiling.
3. Considering the internal spring, it is important to turn the socket pipe in the direction of the arrow and that the fingerscrews (!) are faced forward.
4. Place the socket pipe in the wall bracket and fasten the square-guide with the included screws.
5. The extraction arm can now be turned 360° in the column.

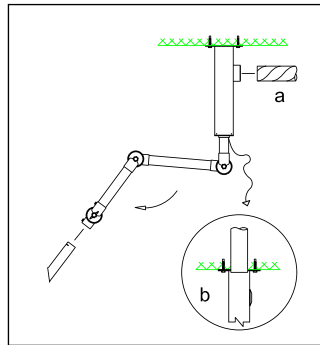**Fig. 2**

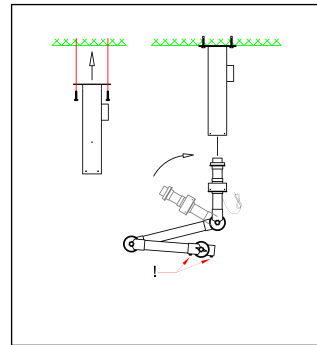1. For side connection (Fig. 2a) the main ventilation duct (80mm dia.) can be connected directly to the outlet on the side of the column.
2. In an ESD area a potential equalizing must be ensured for the surrounding working area. The earthing wire with the 1MΩ resistance must be connected to the central earthing for the furniture and tools. Take care not to break the 1MΩ resistance.
3. Unfold the extraction arm in the direction of the arrow.
4. Place the connection pipe of the accessory in the joint. Turn the small slit in the connection pipe towards the fingerscrew. When screwed in the fingerscrew will hold the accessory in place in the joint.
5. Finally make sure the damper is open.
6. The extraction arm is now ready for use.

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All rights reserved
Alsident® System A/S