HM  
gas

# Alsident® System 100

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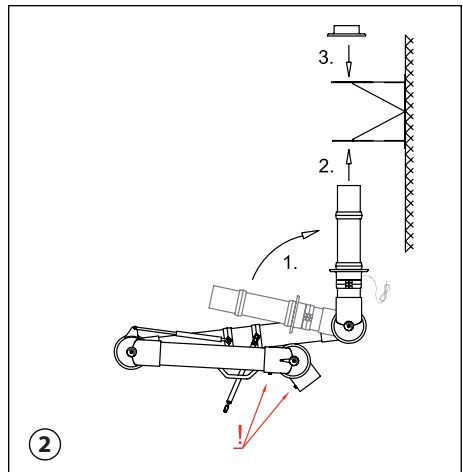
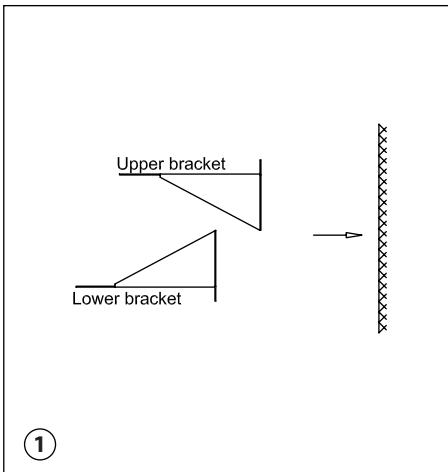
UK

**Fig. 1 (EX/ESD):**

1. The wall bracket is in two parts. Mount the bracket solidly on the wall or column and close together. When mounting the wall bracket be aware of the considerable strain the operation of the arm will transmit to the wall bracket.
2. Please note ! If the wall or column is made of a conductive material the connection to the wall bracket must be non-conductive.
3. In an EX-area the wall bracket must be of stainless steel.

**Fig. 2 (EX/ESD):**

For mounting in an ESD/EX working area the earthing wire on the socket pipe is very important. Take care not to damage it during the mounting.

1. When unfolding the extraction arm, make sure the socket pipe is turned in the direction of the arrow and the finger-screws (!) are faced forward.
2. Place the socket pipe in the bracket and fasten the flange to the bracket with screws.
3. Place the upper flange on the socket pipe and fasten it to the bracket with screws.
4. The two flanges are not fastened to the socket pipe therefore the extraction arm can be turned 190° in the wall bracket.

*NB! Gas springs are considered as wearing parts.*

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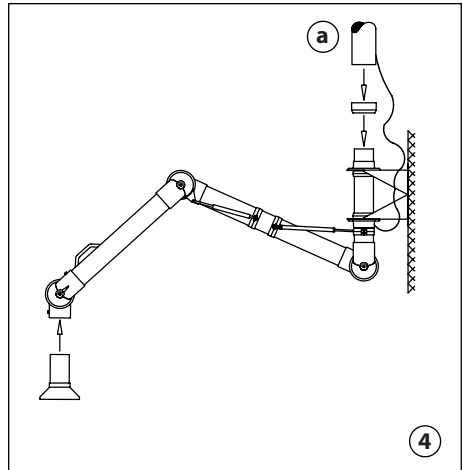
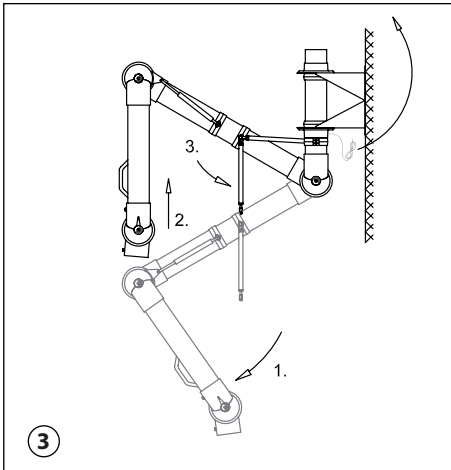
UK

**Fig. 3 (EX):**

1. Unfold the extraction arm in the direction of the arrow (1).
2. Push the extraction arm towards vertical position. This enables the gas spring to be mounted (2).
3. Turn the gas spring up towards the fitting on the socket pipe and fasten it with the fingerscrew (3).
4. In an explosive working area a potential equalizing must be ensured for the ventilation system. It is therefore important to bond all the conductive parts of the system and earthing them.

The extraction arm must be connected to the ventilation duct with the earthing wire pre-attached to the extraction arm. (fig. 4a) Alsident® System A/S recommends the earthing of the extraction arm to be carried out by personnel experienced in EX areas.

**Fig. 4 (EX):**

1. Connect the extraction arm to the main ventilation duct with the Alsident® reducer.
2. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small fingerscrew. When screwed in the fingerscrew will hold the accessory in place in the joint.
3. Finally make sure the damper is open.
4. The extraction arm is now ready for use.

*Approved for use in explosive areas and marked*

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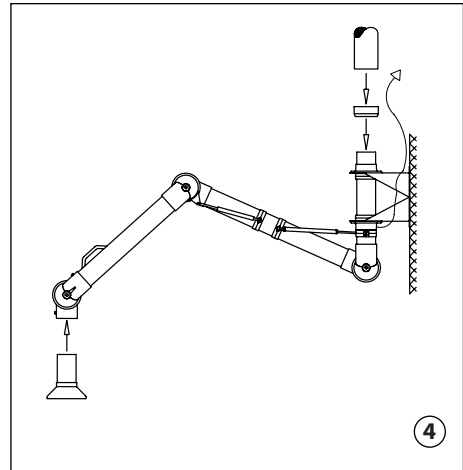
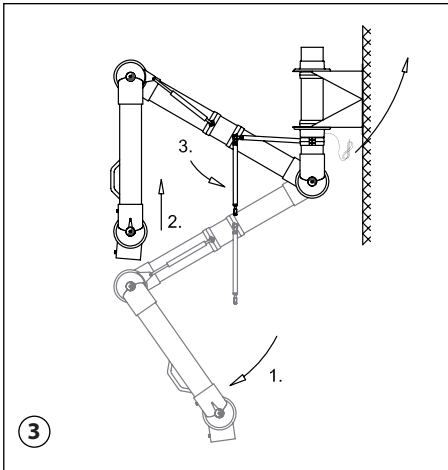
UK

**Fig. 3 (ESD):**

1. Unfold the extraction arm in the direction of the arrow (1).
2. Push the extraction arm towards vertical position. This enables the gas spring to be mounted (2).
3. Turn the gas spring up towards the fitting on the socket pipe and fasten it with the fingerscrew (3).
4. In an ESD area a potential equalizing must be ensured for the surrounding working area. The earthing wire pre-attached to the extraction arm must be replaced with the one in the plastic bag marked »ESD«. The new earthing wire with a 1MΩ resistance must be connected to the central earthing for the furniture and tools (fig. 4). When turning the extraction arm be aware not to break the 1MΩ resistance on the earthing wire. Be aware to break any conductive contact between the bracket and the building/ furniture and main ventilation system.

**Fig. 4 (ESD):**

1. Connect the extraction arm to the main ventilation duct with the Alsident® reducer.
2. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small fingerscrew. When screwed in the fingerscrew will hold the accessory in place in the joint.
3. Finally make sure the damper is open.
4. The extraction arm is now ready for use.

Approved according to standard IEC 61340-5-1:2003.  
For further information, see Technical Information at:  
[www.alsident.com](http://www.alsident.com).

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