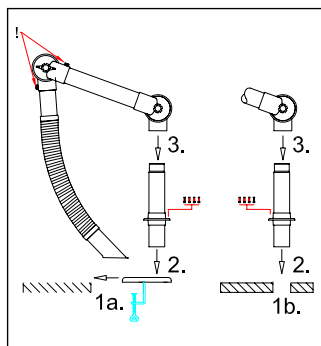
# ALUMINIUM (AL)


## UK – Flex – TM

**Fig. 1**

On account of the packaging the extraction arm is delivered in two parts.

- 1a. Mount the table bracket on the edge of the table and tighten the clamps.
- 1b. Or make a hole in the table with a diameter of 60 mm where the extraction arm is to be placed.
2. Place the socket pipe in the hole of the table/bracket and fasten the flange with the included screws.
3. Place the joint of the extraction arm on the socket pipe. Be aware that the spring goes into the pipe and the fingerscrews (!) face forward.

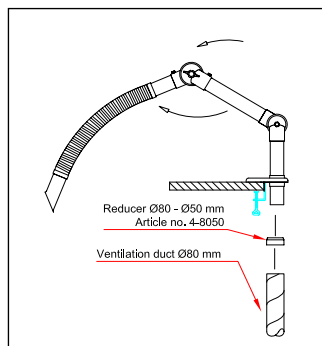
**Fig. 2**

1. Connect the extraction arm to the main ventilation duct with the Alsident® reducer.
2. Unfold the extraction arm in the direction of the arrows.
3. Finally make sure the damper is open.
4. The extraction arm is now ready for use.

- Twist the hose counter-clockwise to increase tension.
- Twist only until the tension required has been reached.
- The hose might break if the tension is too high (when moving the hose).
- Hoses longer than standard length might not be able to stay in position. This also applies if the suction pen is replaced by a hood.
- The hose is a wearing part with limited durability.
- Hoses that have been twisted too much will not be replaced.

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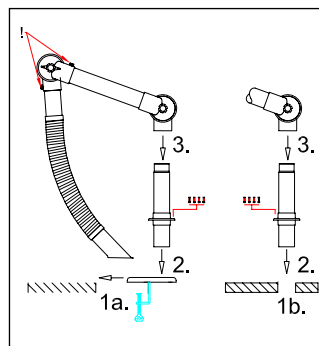
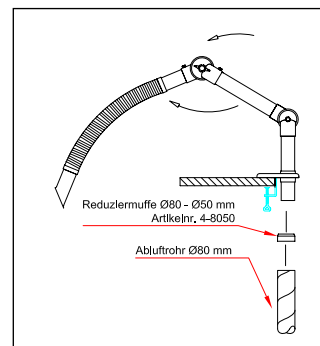
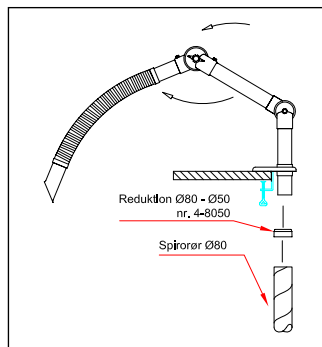
## UK – Flex – TM

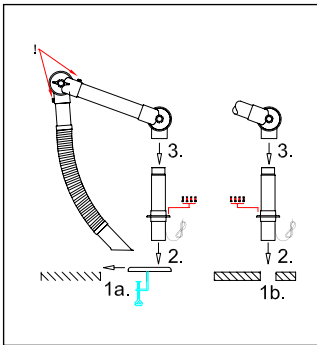
**Fig. 1**

On account of the packaging the extraction arm is delivered in two parts.

For mounting in an ESD working area the earthing wire on the socket pipe is very important. Take care not to damage it during the mounting.

- 1a. Mount the table bracket on the edge of the table and tighten the clamps.
- 1b. Or make a hole in the table with a diameter of 60 mm where the extraction arm is to be placed.
2. If the table is of a conductive material the connection to the extraction arm/bracket must be non-conductive.
3. Place the socket pipe in the hole of the bracket/table and fasten the flange to the bracket/table.
4. Place the joint of the extraction arm on the socket pipe. Be aware that the spring goes into the pipe and the fingerscrews (!) face forward.

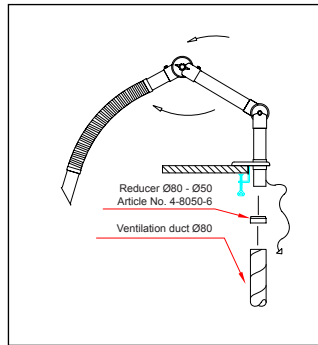
**Fig. 2**

1. Connect the extraction arm to the main ventilation duct with the Alsident® reducer.
2. In an ESD area a potential equalizing must be ensured for the surrounding working area. The earthing wire with the 1MΩ resistance must be connected to the central earthing for the furniture and tools. Take care not to break the 1MΩ resistance.
3. Unfold the extraction arm in the direction of the arrows.
4. Finally make sure the damper is open.
5. The extraction arm is now ready for use.

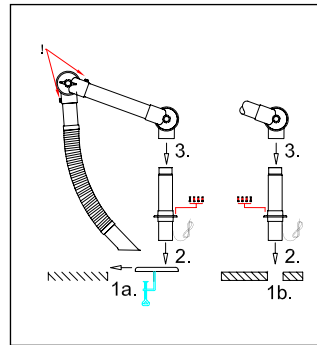- Twist the hose counter-clockwise to increase tension.
- Twist only until the tension required has been reached.
- The hose might break if the tension is too high (when moving the hose).
- Hoses longer than standard length might not be able to stay in position. This also applies if the suction pen is replaced by a hood.
- The hose is a wearing part with limited durability.
- Hoses that have been twisted too much will not be replaced.

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